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## 7. RÉFÉRENCES

- 7.1 Association Française De Normalisation (AFNOR). Validation Certificate N.:3M-01/2-09/89A. 3M Petrifilm Coliform Test. Application for the Enumeration of Coliform Colonies With and Without Gas (as Compared to VRBL Method).
- 7.2 AFNOR. Validation Certificate N.:3M-01/2-09/89B. 3M Petrifilm Coliform Test. Application for the Enumeration of Coliform Colonies With Gas (as Compared to MPN Method).
- 7.3 AFNOR. Validation Certificate N.:3M-01/2-09/89C. 3M Petrifilm Coliform Test. Application for the Enumeration of Thermotolerant Coliform Colonies With and Without Gas (as Compared to VRBL 44.5°C Method).
- 7.4 American Public Health Association. 1992. Standard Methods For the Examination of Dairy Products, 16th ed. APHA, Washington, DC. pp. 248, 251-252, 261-262.
- 7.5 American Public Health Association. 1992. Compendium of Methods for the Microbiological Examination of Foods, 3rd ed. APHA, Washington, DC. pp. 61, 80-89, 338.
- 7.6 AOAC International. 2000. Official Method 986.33. Official Methods of AOAC International. 17<sup>th</sup> ed. AOAC International. Gaithersburg, MD.
- 7.7 AOAC International. 2000. Official Method 989.10. Official Methods of AOAC International. 17<sup>th</sup> ed. AOAC International. Gaithersburg, MD.
- 7.8 AOAC International. 2000. Official Method 991.14. Official Methods of AOAC International. 17<sup>th</sup> ed. AOAC International. Gaithersburg, MD.
- 7.9 Baylis, C. L., C de W. Blackburn and S. B. Pettit, 1994. Evaluation of Petrifilm Plates for the Enumeration of Total Aerobic Flora, Coliforms, and *E. coli*, in Foods. Leatherhead Food RA..

- 7.10 Curiale, M. S., T. Sons, D. McIver, J. S. McAllister, B. Halsey, D. Roblee and T. L. Fox. 1991. Dry Rehydratable Film for Enumeration of Total Coliforms and *Escherichia coli* in Foods: Collaborative Study. J. AOAC. **74**:635-648.
- 7.11 Curiale, M. S., P. Fahey, T. L. Fox and J. S. McAllister. 1989. Dry Rehydratable Films for Enumeration of Coliforms and Aerobic Bacteria in Dairy Products: Collaborative Study. J. AOAC. **72**:312-318.
- 7.12 Food and Drug Administration. 1998. Bacteriological Analytical Manual, 8th ed. Rev. A.. AOAC International, Gaithersburg, MD. 4.11, Appendix 1.10 and Appendix 3.64 (R11).
- 7.13 Ginn, R. E., V. S., Packard and R. L. Fox. 1986. Enumeration of Total Bacteria and Coliforms in Milk by Dry Rehydratable Film Methods: Collaborative Study. J. AOAC. **69**:527-531.
- 7.14 International Dairy Federation (IDF). 1996. International IDF Standard 122C:1996. Milk and Milk Products - Preparation of Samples and Dilutions for Microbiological Examination. Brussels, Belgium.
- 7.15 Matushek, M. G., M. S. Curiale, J. S. McAllister and T. L. Fox. 1992. Comparison of Various Plating Procedures for the Detection and Enumeration of Coliforms in Ice Cream and Ice Milk. J. Food Protection. **55**:113-115.
- 7.16 U.S. Department of Health and Human Services. 1999. Garde "A" Pasteurized Milk Ordinance: Recommendations of the Public Health Service. Publication No. 229. Washington, DC. US Gov't Print Office. 1999 Revision.